

pH METER, pH 3310, WTW

1. Open the box (Fig. 1).
2. Connect the electrode according the picture (Fig. 2).
3. Take out pH meter and pH electrode SenTix 41 (allowed to measure also temperature) (Fig. 3).
4. Place the electrode to the holder (the electrode could be hold in the hand too).
5. Pull off the electrode cover contained 3 M KCl.
6. Wash the electrode by DEMI water.
Switch on the pH meter and immerse the electrode to the measured solution.
7. On the display you see the pH value and the liquid temperature.
In the end of measurement wash the electrode with DEMI water. keep it in 3 M KCl.
8. Switch off the instrument.



Fig. 1: Box with pH Meter.



Fig. 2: Connection of pH electrode.



Fig. 3: Connected pH Meter.

NOTE: If you are not experienced in measuring pH or if you need any assistance or advice, contact Kateřina Tmejová using email: Katerina.Tmejova@ceitec.vutbr.cz