

ANALYTICAL BALANCE ABT 220-5DNM, KERN

1. Move the plastic cover from the balance.
2. Switch-On the balance by quick pushing of the button ON/OFF.
3. There should be a value zero on display.
4. Open glass doors on both sides to be able to work with two hands.
5. Place the plastic bowl/chemical glass. Close the glass doors and push the button TARE.
6. Carefully add the powder to the plastic bowl by the plastic spoon or micro spatula (picture 1a, b).




Picture 1a – Position of the plastic bowl




Picture 1b – Weighing

7. If you want to subtract the value on balance, always close the glass doors and wait few seconds until the value is settled.
8. To remove the plastic bowl, use the tweezers (picture 2).
9. If you make some disorder, carefully brush the mess out on a paper (picture 3, 4).
10. After you finish the weighing, push the OFF button.
11. Put back the plastic cover.

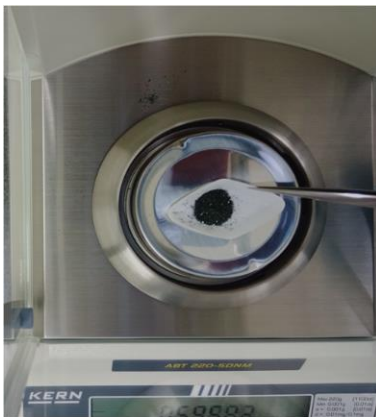

Picture 2 – Removing by tweezers

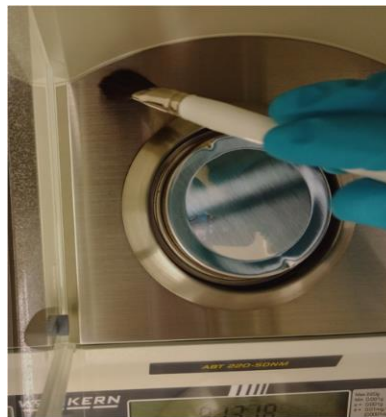

Picture 3 – Cleaning



Picture 4 – Equipment for balance